

Topic

This Technical Note details the new location of Color Tables, Automator Menu Items and Automator Subroutines in Openlab 5 and describes how to move older automations and color tables to the new location.

Discussion

In Openlab 5 the “Color Tables”, “Automator Menu Items” and “Automator Subroutines” folders are now stored in “~/Library/Application Support/Openlab/” (see Fig. 2). Previous version of Openlab stored these files in the application folder (see Fig. 1 below). The new location makes it much easier to maintain these folders when upgrading Openlab. When installing new versions of Openlab it will no longer be necessary to copy automations and color tables into the new Openlab application folder.

To get new or existing automations or color tables to appear in Openlab 5 move or save the files into the appropriate folder located in “~/Library/Application Support/Openlab/” (see Fig. 2).

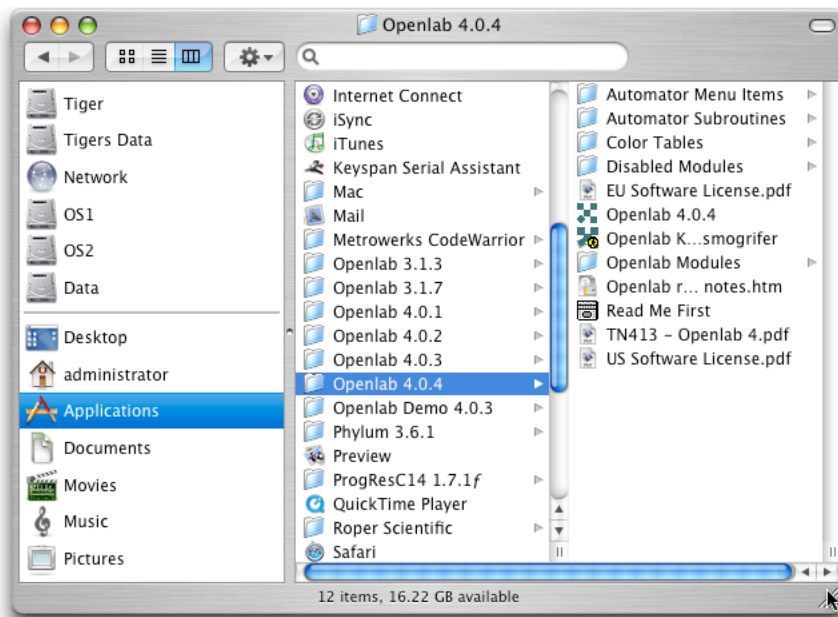


Figure 1. The old location of Color Tables, Automator Menu Items, and Automator Subroutines

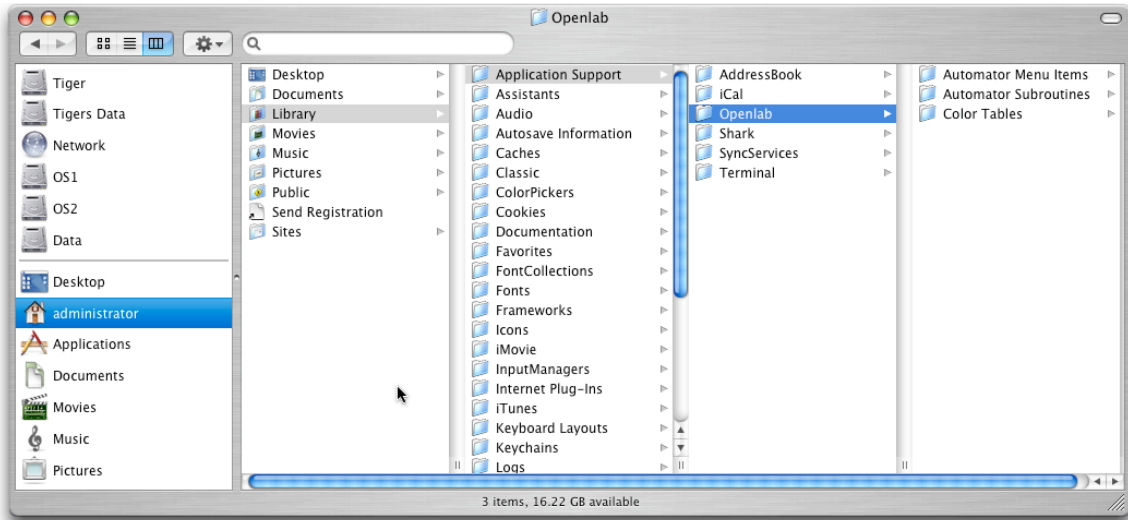


Figure 2. New location of the folders in Application Support. Move old automations and color tables to these folders to load them in Openlab 5.